

Motion Capture Marker Set

Product Description

Motion Capture Marker Set Product Description

v 1.1

October 2021

All rights reserved.

Copyright © Tobii Pro AB (publ)

The information contained in this document is proprietary to Tobii Pro AB. Any reproduction in part or whole without prior written authorization by Tobii Pro AB is prohibited.

Products that are referred to in this document may be either trademarks and/or registered trademarks of the respective owners. The publisher and the author make no claim to these trademarks.

While every precaution has been taken in the preparation of this document, the publisher and the author assume no responsibility for errors or omissions, or for damages resulting from the use of information contained in this document or from the use of programs and source code that may accompany it. In no event shall the publisher and the author be liable for any loss of profit or any other commercial damage caused or alleged to have been caused directly or indirectly by this document.

Content subject to change without notice.

Please visit [Tobii Pro Connect](#) for updated versions of this document.

Table of Contents

1 Overview	4
1.1 Motion Capture Marker Sets	4
1.1.1 Marker Set #1	5
1.1.2 Marker set #2	5
2 Motion capture coordinates	6
2.1 Marker Set #1	6
2.2 Marker Set #2	6
3 Safety	7
3.1 California Proposition 65 warning	7

1 Overview



The Motion Capture Marker set makes it easy to integrate eye tracking data from Tobii Pro Glasses 3 with a motion capture system. For additional information about supported systems, visit our [Motion Capture product page](#).


The set consists of marker holders that attach to the side of the Pro Glasses 3 frame. There are three reflective markers on each holder. The motion capture cameras track the movement of these markers as the subject moves around wearing the Pro Glasses 3. This allows you to stream data from both systems into the motion capture software in real time. The eye tracking and gaze vector data can then be visualized and exported to other analysis software, such as MATLAB.



You can track up to four Pro Glasses 3 at the same time by using multiple marker sets.

1.1 Motion Capture Marker Sets

There are two Marker Sets available, each has different geometry between their reflective markers.

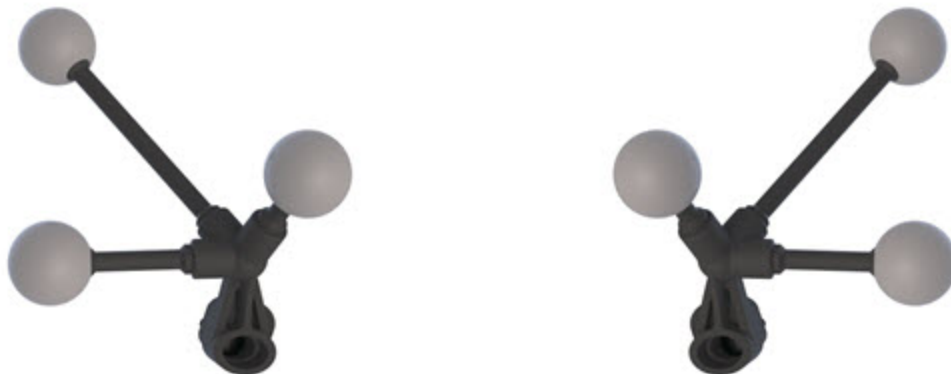

1.1.1 Marker Set #1

Motion Capture Marker Set #1 contains two holders that mount onto Pro Glasses 3. Each holder consists of three passive, reflective markers with a fixed geometry.

1.1.2 Marker set #2

Motion Capture Marker Set #2 enables tracking of two Pro Glasses 3 at the same time when combined with Marker Set #1. The set holds the markers in a different geometry than Marker Set #1. This makes it possible for the motion capture system to distinguish between the different Pro Glasses 3 being tracked.



By combining marker sets, up to four Pro Glasses 3 can be tracked in the same motion capture system.

2 Motion capture coordinates

The coordinates for the reflective markers are listed in the tables below.

2.1 Marker Set #1



	X	Y	Z		X	Y	Z
R1	65.08mm	29.44mm	-1.86mm	L1	-65.08mm	29.44mm	-1.86mm
R2	82.92mm	29.44mm	-52.76mm	L2	-82.92mm	29.44mm	-52.76mm
R3	96.54mm	10.56mm	-18.65mm	L3	-96.54mm	10.56mm	-18.65mm

2.2 Marker Set #2



	X	Y	Z		X	Y	Z
R1	58.25mm	56.40mm	9.39mm	L1	-58.25mm	56.40mm	9.39mm
R2	77.39mm	22.49mm	-34.84mm	L2	-77.39mm	22.49mm	-34.84mm
R3	87.14mm	11.22mm	-15.29mm	L3	-87.14mm	11.22mm	-15.29mm

3 Safety

3.1 California Proposition 65 warning



WARNING: This product can expose you to Carbon Black, which is known to the State of California to cause cancer. For more information, visit [P65Warnings.ca.gov](https://www.p65warnings.ca.gov).



Copyright ©Tobii AB (publ). Not all products and services offered in each local market. Specifications are subject to change without prior notice. All trademarks are the property of their respective owners.

Support for Your Tobii Pro Device

Get Help Online

Visit Tobii Pro Connect for help with your Tobii Pro device. It contains the latest information about contacting Support, links to our Learning Center, and much more. Visit connect.tobii.com.

Contact Your Solution Consultant or Reseller

For questions or problems with your product, contact your Tobii Pro sales representative or authorized reseller for assistance. They are most familiar with your personal setup and can best help you with tips and product training.

Visit tobii.com/contact.